

## Picrosirius Red Staining Of Cardiac Muscle Following Phosphomolybdic Acid Treatment

This method eliminates cytoplasmic staining, revealing thin septa and collagen fibers clearly.

### Fixation

Bouins' Fixative

### Sections

Paraffin @ 7 microns.

### Staining

1. Deparaffinize and hydrate to distilled water.
2. Wash in running tap water, 10 minutes.
3. Rinse in distilled water.
4. Treat in Phosphomolybdic Acid 0.2% Aqueous (1-5 minutes)
5. Stain in Sirius Red, 0.1% in Saturated Picric Acid , for 90 minutes.
6. Wash for 2 minutes in 0.01 N Hydrochloric Acid .
7. Rinse in 70% Alcohol for 45 seconds.
8. Dehydrate, clear and mount in Permount

### Results

Septa          Red

#### 订购信息

H26357	天狼猩红溶液（0.1%，溶于饱和苦味酸）	100ml
--------	----------------------	-------